



1 Detection using VIDAS LSX

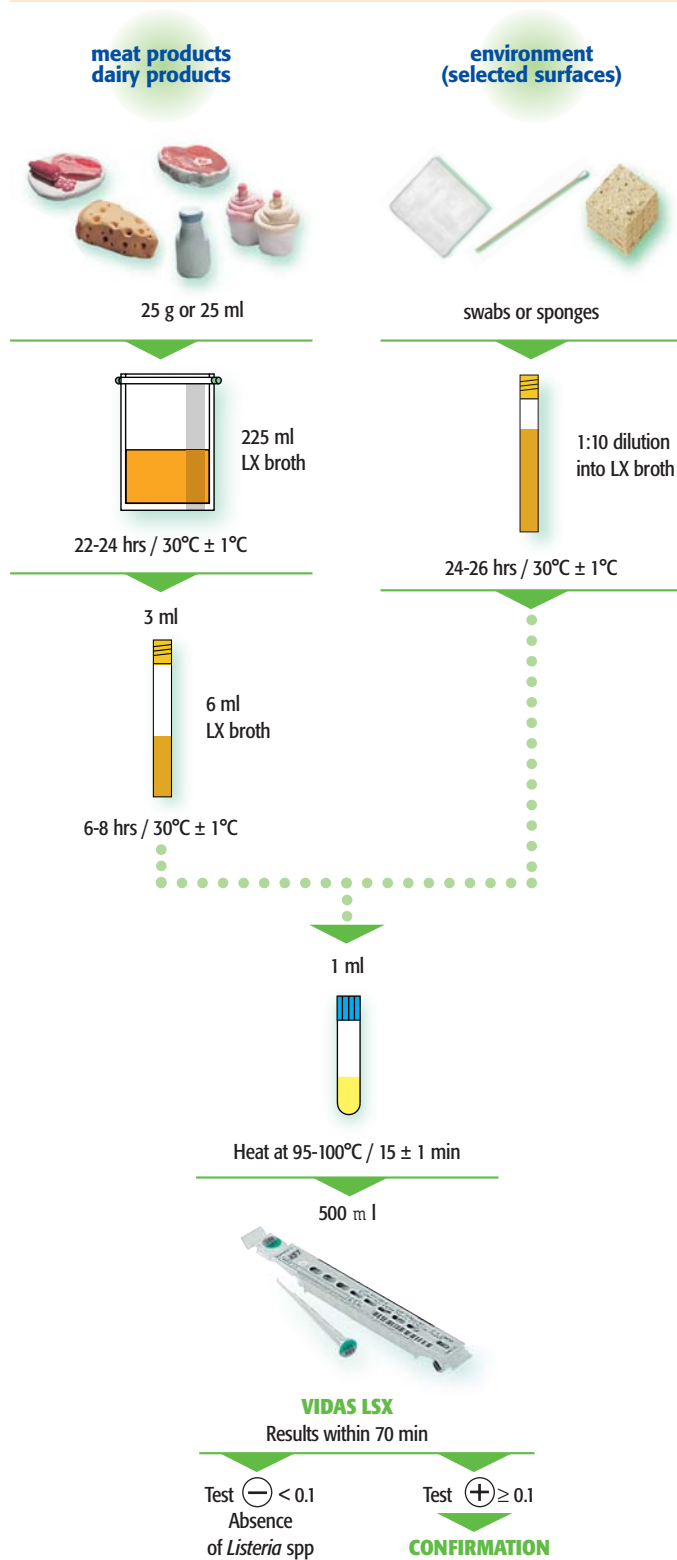

2 Confirmation

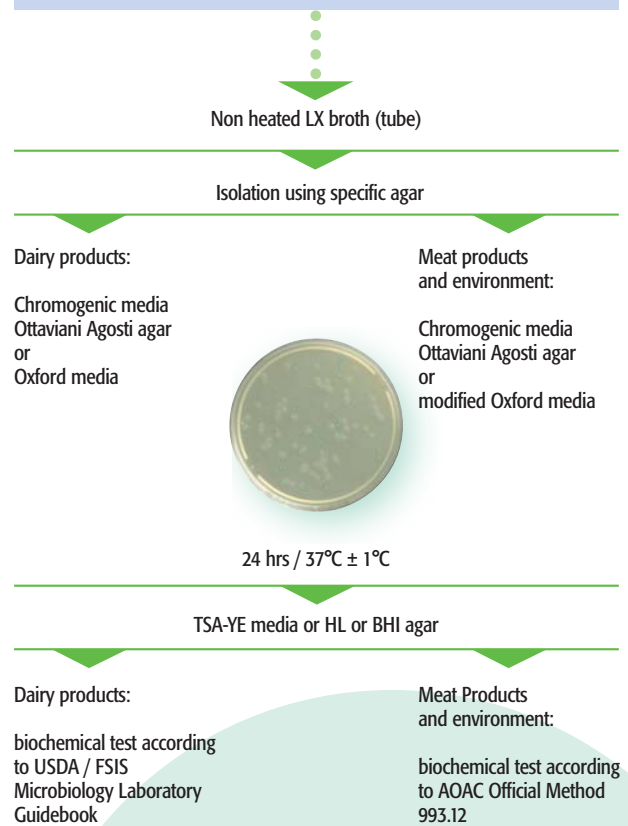

Listeria Spp Xpress range

42626	Listeria Xpress broth	6 x 225 ml
42635	Listeria Xpress broth	6 x 100 ml
42628	Listeria Xpress broth	3 x 3 l
42117	Listeria Xpress broth	20 tubes / 6 ml
42120	Listeria Xpress broth	20 tubes / 10 ml
43560	Oxford agar	20 plates
43641	Ottaviani Agosti agar (OAA)	20 plates
10300	API® Listeria	10 strips
30224	VIDAS® LSX	60 tests


For further details, consult the product package insert.

VIDAS LIS


Detection and confirmation of *Listeria* spp

Protocol validated by AFNOR n° BIO-12/2-06/94 for all food products and environment

1 detection using VIDAS LIS

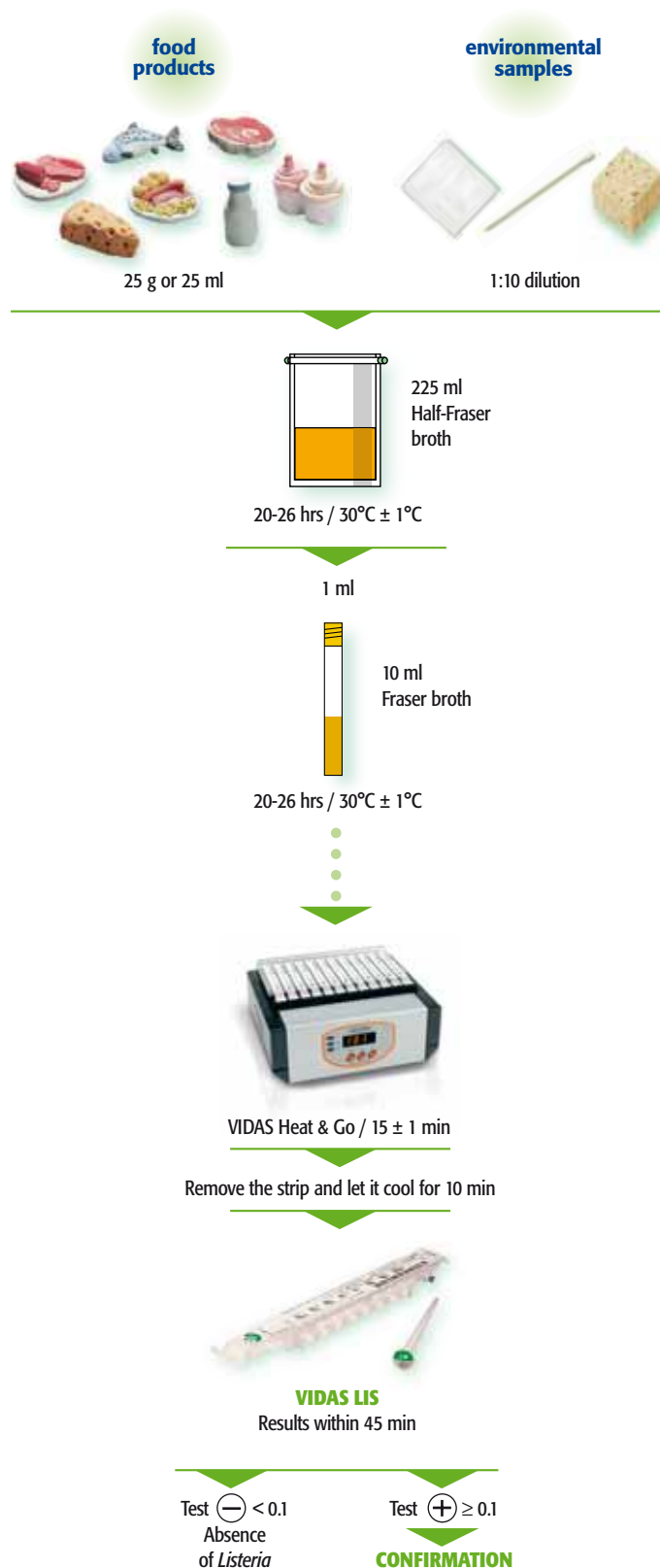

2 confirmation with reference method

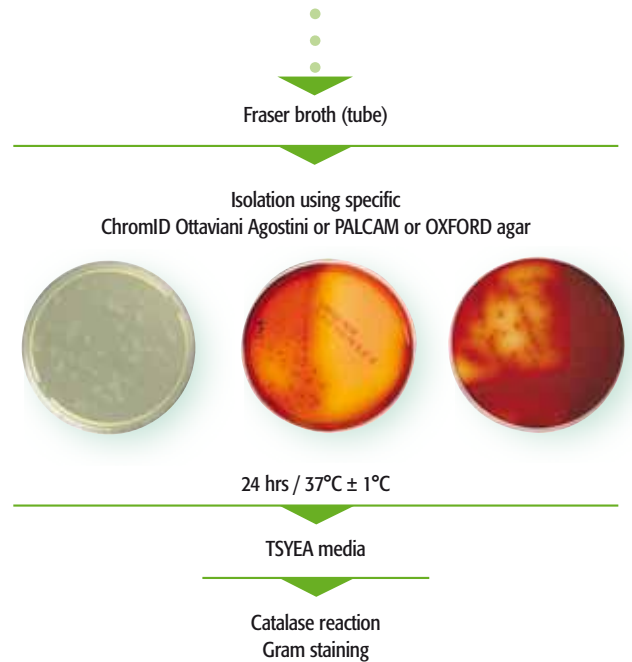

Listeria range

42615	Half-Fraser broth (Bottle 300 ml)	4 x 225 ml
42048	Half-Fraser broth (Bottle 250 ml)	6 x 225 ml
42112	Half-Fraser broth	20 tubes 10 ml
42627	Half-Fraser broth	3 x 3 l
42727	Half-Fraser broth (Minibags)	10 x 225 ml
42072	Fraser broth	20 tubes 10 ml
43559	Palcam agar	20 plates
43560	Oxford agar	20 plates
43641	ChromID Ottaviani Agostini Agar	20 plates
30700	VIDAS Listeria (LIS)	60 tests

For further details, consult the product package insert.



bioMérieux S.A.
69280 Marcy l'Etoile
France
Tel. (33) 04 78 87 20 00
Fax (33) 04 78 87 20 90

www.biomerieux.com
www.biomerieux-industry.com

VIDAS[®] LSX

Detection and confirmation of *Listeria* spp

VIDAS NEXT DAY


Protocol validated by AFNOR n° BIO - 12/12-07/04 for selected food

1 detection using VIDAS LSX

meat products
dairy products
vegetables



25 g or 25 ml

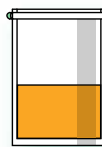

225 ml
LX broth

22-24 hrs / 30°C ± 1°C

3 ml



6 ml
LX broth

6-24 hrs / 30°C ± 1°C

1 ml



Heat at 100°C / 15 ± 1 min

500 µl



VIDAS LSX

Results within 70 min

Test \ominus < 0.1
Absence
of *Listeria* spp

Test \oplus \geq 0.1

CONFIRMATION

2 confirmation

Non heated LX broth (tube)

Isolation using specific agar
Chromogenic media or Palcam agar or Ottaviani Agosti agar or Oxford agar



24 hrs / 37°C ± 1°C

TSYEA media

Catalase reaction
Gram staining

Listeria Spp Xpress range

42626	Listeria Xpress broth	6 x 225 ml
42628	Listeria Xpress broth	3 x 3 l
42117	Listeria Xpress broth	20 tubes
43559	Palcam agar	20 plates
43560	Oxford agar	20 plates
43641	Ottaviani Agosti agar (OAA)	20 plates
10300	API [®] Listeria	10 strips
30224	VIDAS LSX	60 tests

For further details, consult the product package insert.